产品名称 Product Name:	CHO CD Medium 01-BA1201
货号 Cat. No.	BA1201
规格 Amount:	10L、100L、500L、定制 Customized
形式 Formulation:	干粉 Powder
储存温度 Storage:	2 ~ 8°C
有效期 Validity period:	24个月 (失效日期见产品包装) 24 months (Expiration date on product packaging)

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Description

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CHO CD Medium 01-BA1201 provides a rich environment for high density CHO cell growth. CHO CD Medium 01-BA1201 is a serum-free, animal component-free, chemically defined medium developed for CHO suspension culture for expression of antibodies and protein products. CHO suspension cultures can be subculture directly into CHO CD Medium 01-BA1201 from serum-supplemented or serum-free medium with little or no adaptation. CHO CD Medium 01-BA1201 is formulated without hypoxanthine, thymidine and L-glutamine, suitable for DHFR, glutamine synthetase (GS System) screening system.

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Use aseptic technique when handling or supplementing this medium. This product is for research or for further

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Caution: Not for human or animal therapeutic use. Uses other than the intended use may be a violation of local law.

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Read the Material Safety Data Sheets (MSDS) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

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Add 80~90% of the final volume of water for injection to a suitable, clean container, and adjust the water temperature to 25~35°C;

2. 缓慢加入 21.2g /L 的 CHO CD Medium 01-BA1201 干粉培养基,缓慢加入水中,充分搅拌 10 分钟,溶液不澄清; Slowly add 21.2 g/L CHO CD Medium 01-BA1201 powder, stir and mix for 10 minutes, the solution does not clear;

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Slowly add 5 mol/L NaOH adjust the pH to 9.05 ± 0.1 and stir 30 minutes;

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Slowly add 6 mol/L HCl adjust the pH to 7.00 ± 0.1 and stir 10minutes, add 3.575 g/L NaHCO₃ and stir 10 minutes until dissolved;

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Slowly add 6 mol/L HCl to adjust the pH to 7.0 ± 0.1 and stir 10 minutes;

6. 加注射用水定容到最终的配液体积，继续搅拌 1-2 分钟。
Add water for injection to the desired final volume, stir 1-2 minutes;

7. 通过 0.22 μ m 孔径滤膜过滤除菌，2-8°C 避光保存。
Sterilize by filtration using pore size 0.22 μ m filter, 2 ~ 8°C protect from light.

- CHO CD Medium 01-BA1201 培养基的使用需要无菌；
Use CHO CD Medium 01-BA1201 under sterile condition;
- 产品不含 L-谷氨酰胺；使用前按需添加 L-谷氨酰胺；
This product is formulated without L-glutamine, add L-glutamine as needed before use;
- 不推荐使用抗生素；
Antibiotics are not recommended;
- 在 2 ~ 8°C 避光干燥保存。
2~8°C, keep dry and protect from light.

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Transfer all the cell suspension in the centrifuge tube to a 125 mL shake flask containing 15 mL of preheated complete CHO CD Medium 01-BA1201 medium and dilute to the required cell density;

4. 在含有 5 ~ 8% CO₂，37°C，加湿的培养箱或摇床进行培养，培养时拧松瓶盖或使用通气盖以进行气体交换；
Culture in incubator contains 5% ~ 8% CO₂, 37°C humidified air; Unscrew the cap or use the aeration cap for gas exchange during incubation.

5. 细胞复苏后培养 2 ~ 5 天处于对数生长期中期时传代。在进行其它实验之前，复苏的细胞至少应进行三次传代。
Cells should be cultured for 2 to 5 days after cell recovery and passed on in the middle stage of logarithmic growth. The cells should be passed on at least three times before other experiments.

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Related Products

货号 Cat.No.	品名 Product Name
BA1202	CHO CD Feed 01-BA1202
BA1203	CHO CD Feed 02-BA1203
BA1204	CHO CD Feed 01 Plus-BA1204

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Explanation of Symbols and Warnings

		
供 GMP 制造 For GMP Manufacturing	有效期至 Expiration Date	储存温度 Store Temperature
		
批号 Lot Number	干燥保存 Keep dry	避光保存 Protect from light